



**EDEN LAB**  
LABORATORY SUPPLIES



Authorised Distributor  
For

**RICHTER**  **PTICA**

[contact@edenlab.com.au](mailto:contact@edenlab.com.au)  
[www.edenlab.com.au](http://www.edenlab.com.au)

# UX1 Microscopes


---





# UX1 Microscope Features

- 1** Ergonomically inclined head with eyetubes at 30° places user in the best working position, creating less tension and fatigue. Interpupillary distance is adjustable from 50~75mm.
- 2** Trinocular port with c-mount adapter for camera or tablet connection.
- 3** Double layer mechanical stage 140mm x 132mm with 50mm x 76mm movement range.
- 4** Abbe condenser is pre-centered and non-removable to maintain factory preset. Image contrast is improved with iris diaphragm and filter holder.
- 5** Coaxial coarse & fine focus adjustment on both sides of microscope frame. Coarse tension is adjustable. Upper limit avoids impact between objective and slide.
- 6** LED illumination is energy efficient with long-lasting bulb. Rechargeable batteries allow cordless or corded microscope operation.
- 7** Select from Achromat or Plan Achromat objective lenses.
- 8** Rheostat control for light intensity control.

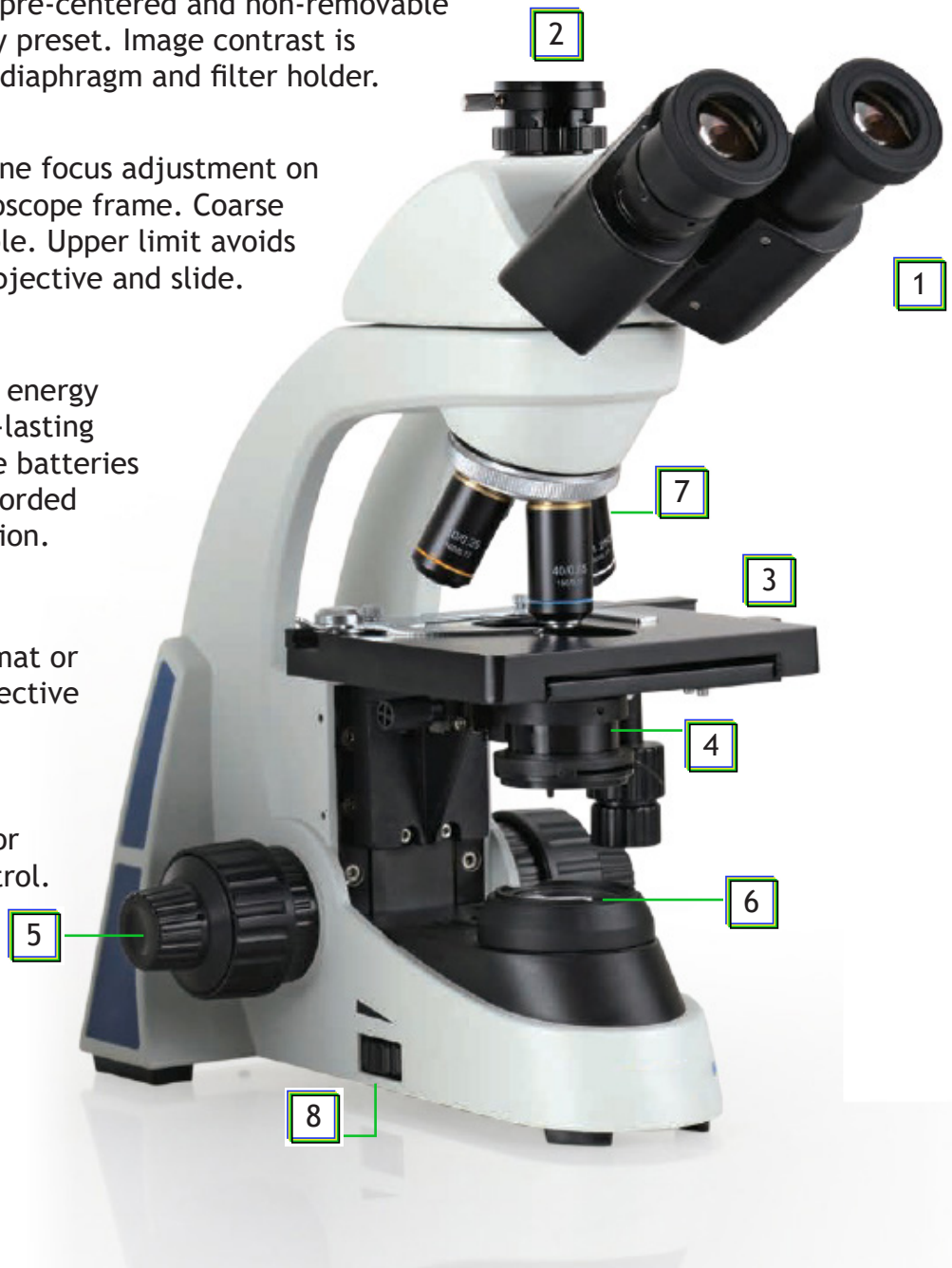



# UX1 Objective Lenses



## Achromat Objectives

Objective	N.A.
4x	0.10
10x	0.25
40xs	0.65
100xs oil	1.25



## Plan Achromat Objectives



Objective	N.A.	W.D.
4x	0.10	15.1mm
10x	0.25	1.5mm
40xs	0.65	0.38mm
100xs oil	1.25	0.09mm





# UX1 Microscope Models



**UX-1B Binocular**

Binocular eyepieces have an interpupillary distance adjustable from 50~75mm.



**UX-1T Trinocular**

Trinocular port for camera attachment.  
Beam splitter ratio = 80:20.



**UX-1D Digital**

3mp or 5mp USB Camera with Software  
for image capture and measurement.





**UX1-LCD Tablet**

8" LCD Touch-Screen Tablet for image  
display and capture.

# UX1 Microscope Specifications



Optical System	Finite Color Corrected Optical System
Viewing Head	30° inclined binocular head. Interpupillary distance adjustable 50-75mm. Left eyetube is +/-5 diopter adjustable.
	30° inclined trinocular head. Interpupillary distance adjustable 50-75mm. Left eyetube is +/-5 diopter adjustable. Beam splitter ratio = 80:20 when engaged.
Eyepieces	Widefield plan eyepiece PL10x, FN18. Retaining ring available for 19mm diameter reticle.
Objective Lenses	Achromat 4x, 10x, 40xs, 100xs oil.
	Plan Achromat 4x, 10x, 40xs, 100xs oil.
Nosepiece	Reversed quadruple nosepiece.
Focus Adjustment	Coaxial coarse and fine focus adjustment with coarse adjustment stop and tension adjustment. Range: 25mm, precision 0.002mm.
Stage	132mm x 140mm double layer mechanical stage with moving range 50mm x 76mm precision 0.1mm.
Condenser	N.A. 1.25 Abbe condenser with iris diaphragm and filter holder.
Illumination	3W LED illuminator with rheostat control, 100V~240V. Includes four AA rechargeable batteries and recharger. Operates plugged in or cordless.
C-Mount Adapters	Focusing 0.35x and 0.5x.

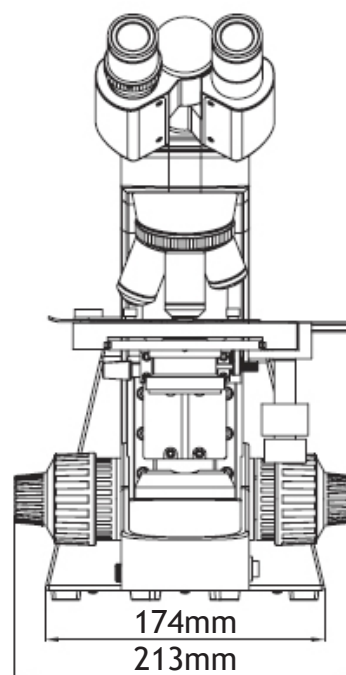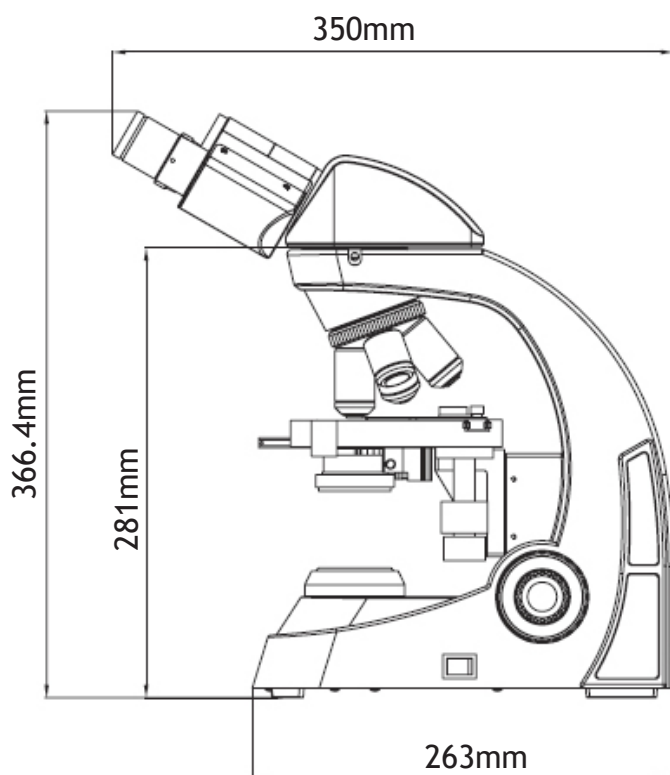

info@richter-optica.com  
www.richter-optica.com

- Autoclaves
- Balances & Scales
- Bath & Circulators
- Centrifuges
- Chillers/ Freeze Dryers
- Desiccators
- Filtration Systems
- Furnace/ Burners
- General Lab Equipment
- Glovebox
- Heating & Cooling blocks
- Heating mantle/ Temperature controls
- Homogenizers
- Hotplate & Stirrers
- Incubators
- Lab Cabinets & Storage
- Laboratory consumables
- Measurement & Testing  
(pH/ DO/ TDS/ Refractometer/Conductivity)
- Microscope (Research)
  - o Life science Microscope
  - o Geology Microscope
  - o Industrial Inspection Microscope
  - o Material Science Microscope
  - o Microscope Cameras
- Microscope (Educational/ Teaching)
  - o Highschool
  - o College
  - o University
- Ovens
- Pipette
- Plasticwares
- Rangehoods/ Clean Bench, BIO Safety Cabinet
- Recirculating Coolers
- Refrigerator/ Freezers
- Rotary Evaporator
- Rotators
- Shakers/ Mixer/ Roller
- Temperature/ Environment Chambers
- Ultrasonic cleaner
- UV Vis Spectrophotometer
- Vacuum Pump

